

## DISCLAIMER

This software has not undergone extensive formal testing at this time. It will be included in the next version of the Analyst® software, at which time it will have been subjected to such testing. SCIEX assumes no liability or responsibility for any loss howsoever caused in connection with using this software.

# Analyst® Software

Version 1.6.3

## Analyst® 1.6.3 Patch for Checksum Issue

This patch fixes the defect where, intermittently, data for one of the samples within the batch is not recorded and that data file might also have an invalid checksum. (AN-1250)

### Where to Get Help

#### Support

- <https://sciex.com/request-support>

### Install the Patch

1. Deactivate the hardware profile and then close the Analyst® software.
2. Double-click the **Analyst 1.6.3 Patch for Checksum Issue.exe** file and then follow the on-screen instructions.

#### Updated Files

The patch updates the Analyst® software file, FMWIFFCompDocNTDriver.dll, in the following folder:

Windows 7, 64-bit operating systems: C:\Program Files (x86)\Analyst\bin

Windows 7, 32-bit operating systems: C:\Program Files\Analyst\bin

## Remove the Patch

1. Open the Control Panel and then click **Programs and Features**.
2. Select **Analyst 1.6.3 Patch for Checksum Issue** from the list and then click **Uninstall**.

## Known Issues and Limitations

None

## Notes on Use

None

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate those manufacturers' products as supplied by SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products, and are SCIEX's sole and exclusive representations, warranties, and obligations. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser, or for any adverse circumstances arising therefrom.

**For research use only.** Not for use in diagnostic procedures.

AB Sciex is doing business as SCIEX.

The trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners.

AB SCIEX™ is being used under license.

© 2018 AB SCIEX



AB Sciex Pte. Ltd.  
Blk 33, #04-06  
Marsiling Ind Estate Road 3  
Woodlands Central Indus. Estate  
SINGAPORE 739256

