

DISCLAIMER

This software has not undergone extensive formal testing at this time. It will be included in the next version of the Analyst® Software, at which time it will have been subjected to such testing. SCIEX assumes no liability or responsibility for any loss howsoever caused in connection with using this software.

Analyst® Software, Version 1.6.3

Analyst® 1.6.3 Patch for Shimadzu LC30 Plate Layout (Rev.2)


The Analyst® 1.6.3 Patch for Shimadzu LC30 Plate Layout enables support for a new plate layout for the Shimadzu SIL-30ACMP and the Shimadzu SIL-30AC autosamplers (JIRA AN-1571):

- 96 deep well plate (Alpha Deep Well MTP 96 for the Shimadzu SIL-30ACMP autosampler and for the Shimadzu SIL-30AC autosampler).

The following are applicable when this new plate is selected for the Shimadzu SIL-30ACMP autosampler or for the Shimadzu SIL-30AC autosampler:

- The locations in the Batch Editor are assigned numerical values, arranged horizontally.
- The Batch Editor supports the “fill down” feature.
- The Batch Editor can export to txt and csv files.
- The Batch Editor can import from txt and csv.

Figure 1: Batch Editor – Locations Tab

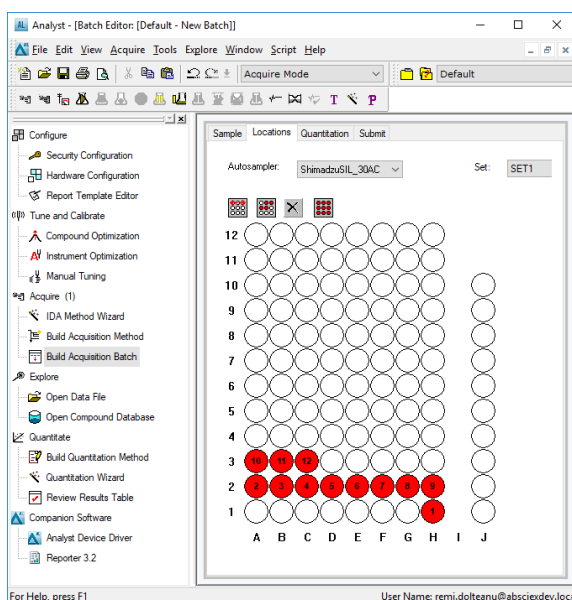

Release Notes

Where to Get Help

Support

- <https://sciex.com/request-support>
- www.sciex.com

Install the Patch

1. Deactivate the hardware profile and then close the Analyst® Software.
2. Uninstall the **Analyst 1.6.3 Patch for Shimadzu LC30 Plate Layout**. Refer to steps 2 through 4 in [Remove the Patch](#).
3. Double-click **Analyst 1.6.3 Patch for Shimadzu LC30 Plate Layout.exe** and then follow the on-screen instructions.

Updated and Added Files

On a computer configured with the Windows 7, 32-bit operating system, the *<path>* refers to <drive:>\Program Files.

On a computer configured with the Windows 7, 64-bit or the Windows 10 operating system, the *<path>* refer to <drive>:\Program Files (x86).

This patch updates the following files in the *<path>*\Analyst\bin folder:

- DDIShimadzu.dll

This patch adds the following file in the *<path>*\Analyst\bin folder:

- AutosamplerDB_SIL30AC_SIL30ACMP_AlphaMTP96.adb

Remove the Patch

1. Deactivate the hardware profile and then close the Analyst® Software.
2. Click **Start > Control Panel**.
3. In the **Large icon** or **Small icons** view, click **Programs and Features**.
4. Select **Analyst 1.6.3 Patch for Shimadzu LC30 Plate Layout** and then click **Uninstall**.

Known Issues and Limitations

None.

Release Notes

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate those manufacturers' products as supplied by SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products and are the sole and exclusive representations, warranties, and obligations of SCIEX. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser or for any adverse circumstances arising therefrom.

GEN-IDV-09-10816-A

For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein are the property of AB Sciex Pte. Ltd., or their respective owners, in the United States and/or certain other countries.

AB SCIEX™ is being used under license.

© 2020 DH Tech. Dev. Pte. Ltd.



AB Sciex Pte. Ltd.
Blk 33, #04-06 Marsiling Industrial Estate Road 3
Woodlands Central Industrial Estate, Singapore 739256