
DISCLAIMER

This software patch has not undergone extensive formal testing at this time. The patch has undergone sufficient testing to prove that it resolves the issue detailed below. This patch will be included in the next version of the Analyst Software, at which time it will have been subjected to complete and formal testing. SCIEX assumes no liability or responsibility for any loss howsoever caused in connection with using this software.

Analyst 1.7.2 Patch for Full User Name (Rev.2)

This patch addresses the following issues:

- Only the last name of the user was shown instead of the full name in the **Full User Name** field of the audit trail. Reports created with a report template also showed only the last name in the **Operator** field (if used). (AN-2221)
- If the full name of the user in the Windows Server Active Directory is longer than 31 characters, then the **Full User Name** shown in the audit trail or in reports could be shown with invalid characters. This issue only occurred in the first release of the Analyst 1.7.2 Patch for Full User Name. (AN-2407)

This patch includes the Analyst 1.7.2 Patch for Screen Lock Issue. If that patch is already installed on the system, then it will be removed during the installation of Analyst 1.7.2 Patch for Full User Name (Rev.2).

The Analyst 1.7.2 Patch for Screen Lock Issue addresses the following issues:

- In Mixed Mode, the dialog to unlock the screen allowed the user to select a domain from a list, and the list was not always complete. Now, the **User Name** field accepts the user name in both Security Account Manager (SAM) and User Principle Name (UPN) formats. (AN-2187)
- In Mixed Mode, users were unable to unlock the screen if their passwords were more than 13 characters long. Passwords up to 20 characters long are now supported. (AN-1194)

Install the Patch

Prerequisites
<ul style="list-style-type: none">• The Analyst 1.7.2 Software is installed.• The previous version of the patch, Analyst 1.7.2 Patch for Full User Name, has been removed if it was already installed. Refer to the section Remove the Patch.

1. Log on to the computer as a user with Administrator privileges.
2. Wait for any acquisitions to complete.
3. Deactivate the hardware profile.
4. Close the Analyst Software.
5. Double-click the **Analyst_1.7.2_Patch_for_Full_User_Name_Rev_2.exe** file.

Note: If the Analyst 1.7.2 Logger for Screen Lock Issue V.2.0 is installed, then it is automatically removed.

6. Follow the on-screen instructions to complete the installation.

Updated Files

The Analyst 1.7.2 Patch for Full User Name (Rev.2) makes the following changes to the Analyst\bin folder.

Note: On 64-bit systems, this folder is in the C:\Program Files (x86)\ folder. On 32-bit systems, it is in the C:\Program Files\ folder.

- AdminConsole.dll (updated)
- Analyst.exe (updated)
- MMSecurity.dll (updated)
- UserManager.dll (updated)
- PEIUtils.dll (updated)

Remove the Patch

1. Log on to the computer as a user with Administrator privileges.

2. Wait for any acquisitions to complete.
3. Deactivate the hardware profile.
4. Close the Analyst Software.
5. Open the Control Panel in Large icons or Small icons mode, and then click **Programs and Features**.
6. Select **Analyst 1.7.2 Patch for Full User Name (Rev.2)** and then click **Uninstall**.
The patch is removed from the program list.

Known Issues and Limitations

- N/A

Release Notes

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate those manufacturers' products as supplied by SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products and are the sole and exclusive representations, warranties, and obligations of SCIEX. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser or for any adverse circumstances arising therefrom.

(GEN-IDV-09-10816-D)

For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd., or their respective owners, in the United States and/or certain other countries (see sciex.com/trademarks).

AB SCIEX™ is being used under license.

© 2021 DH Tech. Dev. Pte. Ltd.



AB Sciex Pte. Ltd.
Blk33, #04-06 Marsiling Industrial Estate Road 3
Woodlands Central Industrial Estate, Singapore 739256