

## DISCLAIMER

This software has not undergone extensive formal testing at this time. It will be included in the next version of the ProMassProcessor Software, at which time it will have been subjected to such testing. SCIEX assumes no liability or responsibility for any loss howsoever caused in connection with using this software.

## ProMassProcessor Software

Version 1.0

## ProMassProcessor 1.0 Patch for Improvements to Inserting Sequence Data

This patch is distributed to users of the ProMassProcessor 1.0 Software. It provides the following improvement:

In version 1.0 of the ProMassProcessor Software, the sequence information was specified on a per sample basis. This patch provides a high-throughput workflow by enabling users to specify the sequence information for multiple samples at a time. The list of sequences can be copied from multiple cells in an Excel spreadsheet and then pasted in the Sequence column in the ProMassProcessor software.

### Where to Get Help

#### Support

<https://sciex.com/request-support>

### Install the Patch

1. Close the ProMassProcessor software.
2. Save **ProMassProcessor 1.0 Patch for Improvements to Inserting Sequence Data.exe** in a folder on the hard drive: <Drive>:\ProMassProcessor 1.0 Patch for Improvements to Inserting Sequence Data\.
3. Navigate to <Drive>:\ProMassProcessor 1.0 Patch for Improvements to Inserting Sequence Data\, double-click **ProMassProcessor 1.0 Patch for Improvements to Inserting Sequence Data.exe**, and then follow the on-screen instructions.

### Updated files

This patch updates the following ProMassProcessor Software files:

- Sciex.ProMassProcessor.exe in the <drive>:\Program Files\SCIEX\ProMassProcessor folder

# Release Notes

---

## Remove the Patch

1. Open Control Panel and then click **Programs and Features**.
2. Select **ProMassProcessor 1.0 Patch for Improvements to Inserting Sequence Data** from the list and then click **Uninstall**.

**Note:** Removal of the ProMassProcessor 1.0 Patch for Improvements to Inserting Sequence Data restores the original ProMassProcessor 1.0 Software.

## Known Issues and Limitations

None

# Release Notes

---

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate those manufacturers' products as supplied by SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products and are SCIEX's sole and exclusive representations, warranties, and obligations. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser or for any adverse circumstances arising therefrom.

For Research Use Only. Not for use in Diagnostic Procedures.

AB Sciex is operating as SCIEX.

The trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners.

AB SCIEX™ is being used under license.

© 2019 AB Sciex



AB Sciex Pte. Ltd.  
Blk 33, #04-06 Marsiling Ind Estate Road 3  
Woodlands Central Indus. Estate SINGAPORE 739256