

## DISCLAIMER

This software has not undergone extensive formal testing at this time. It will be included in the next version of the ProteinPilot™ Software, at which time it will have been subjected to such testing. SCIEX assumes no liability or responsibility for any loss howsoever caused in connection with using this software.

# ProteinPilot™ Software

Version 5.0.2

## ProteinPilot 5.0.2 Patch for UniProt and Issue (Rev.2)

This patch addresses the issue where users are unable to retrieve UniProt annotations. In addition, it also addresses the issue where users are unable to export the Protein Summary and FDR export.

The issue was initially fixed in the release of the ProteinPilot™ 5.0.2 Software. However, UniProt subsequently made additional changes on their website and the UniProt annotations can no longer be retrieved using the ProteinPilot™ Software. UniProt has also changed the xml format that returns an empty sub node, which causes the ProteinPilot™ Software to return an error during the export. This patch fixes both issues.

### Where to Get Help

#### Support

- <http://sciex.com/request-support>

### Install the Patch

**Note:** ProteinPilot™ 5.0.2 Software must be installed on Windows 10, 64-bit systems.

- Double-click **ProteinPilot 5.0.2 Patch for UniProt Issue (Rev. 2).exe** file and then follow the on-screen instructions.

### Updated Files

The patch updates the **progroup.exe**, **ABSciex.ProteinObjects.dll**, **ABSciex.ProteinObjects\_x64.dll** file in the C:\Program Files\AB SCIEX\ProteinPilot folder.

# Release Notes

---

## Remove the Patch

1. Open **Control Panel** and then click **Programs and Features**.
2. Select **ProteinPilot 5.0.2 Patch for UniProt Issue (Rev. 2)** from the list and then click **Uninstall**.

**Note:** Removal of the ProteinPilot 5.0.2 Patch for UniProt Issue (Rev.2) restores the original ProteinPilot™ 5.0.2 Software version.

## Known Issues and Limitations

None

# Release Notes

---

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate those manufacturers' products as supplied by SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products and are the sole and exclusive representations, warranties, and obligations of SCIEX. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser or for any adverse circumstances arising therefrom.

For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein are the property of AB Sciex Pte. Ltd., or their respective owners, in the United States and/or certain other countries. AB SCIEX™ is being used under license.

© 2020 DH Tech. Dev. Pte. Ltd.



AB Sciex Pte. Ltd.  
Blk 33, #04-06 Marsiling Ind Estate Road 3  
Woodlands Central Indus. Estate SINGAPORE 739256