

Release Notes


DISCLAIMER

This software fix has not undergone extensive formal testing at this time. The fix has undergone sufficient testing to prove that it resolves the issue detailed below. This patch will be included in the next version of SCIEX OS, at which time it will have been subjected to complete and formal testing. SCIEX assumes no liability or responsibility for any loss howsoever caused in connection with using this software.

SCIEX OS Version 1.7

SCIEX OS 1.7 Patch for Installation on German OS

This patch is distributed to SCIEX OS customers running SCIEX OS on a computer with Windows configured for German. It addresses the following issue: Customers cannot install SCIEX OS 1.7 for data processing when the Windows OS is configured for German. When the user tries to install the following error is shown. (BLT-1942)

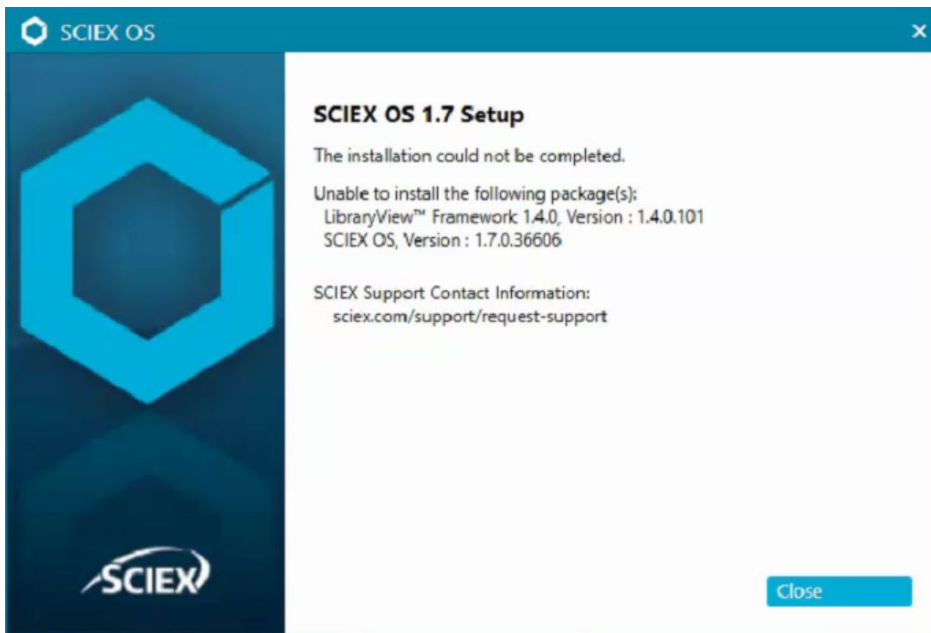

Where to Get Help

Support

- www.sciex.com/support
- www.sciex.com

Release Notes

Install the Patch

1. Save **SCIEX_OS_1.7_Patch_for_Installation_on_German_OS_Issue.zip** in a folder on the hard drive (**D:\SCIEX OS 1.7 Patch for Installation on German OS**) and then extract it.
2. Run **setup.exe** from the SCIEX OS 1.7 package.
Note: Installation of the LibraryView™ Framework fails.
3. Navigate to the folder **D:\SCIEX OS 1.7 Patch for Installation on German OS**, double-click **LibraryView_Framework_x64.msi**, and then follow the instructions.
4. Run **setup.exe** from the SCIEX OS 1.7 package again.
All of the necessary components for SCIEX OS 1.7 are installed.

Updated files

LibraryView™ Framework 1.4.0.124 is installed.

Remove the Patch


N/A

Release Notes


Known Issues and Limitations

(Windows 7) For both English and German systems, the SCIE X OS installation fails if no version of .NET 4.x is installed.

Note: This issue does not occur on computers supplied by SCIE X. (ONYX-8028)

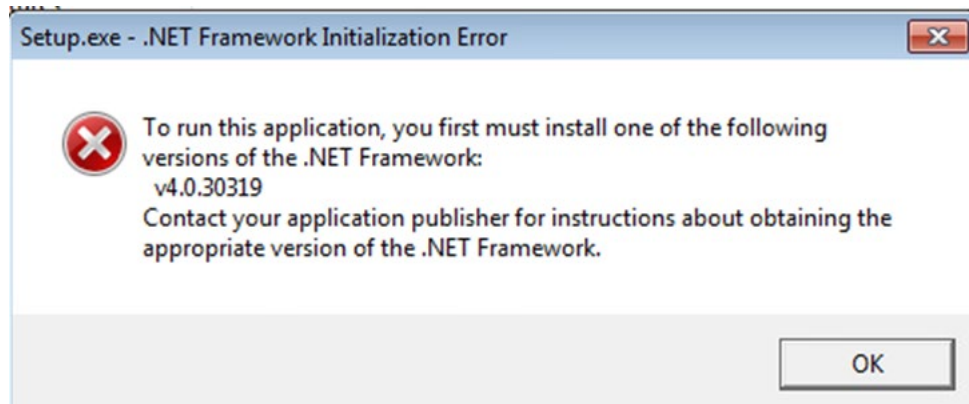

Workaround: If this issue occurs, then run **NDP472-KB4054530-x86-x64-AIIOS-ENU.exe**, located in the SCIE X OS 1.7 **Install** folder.

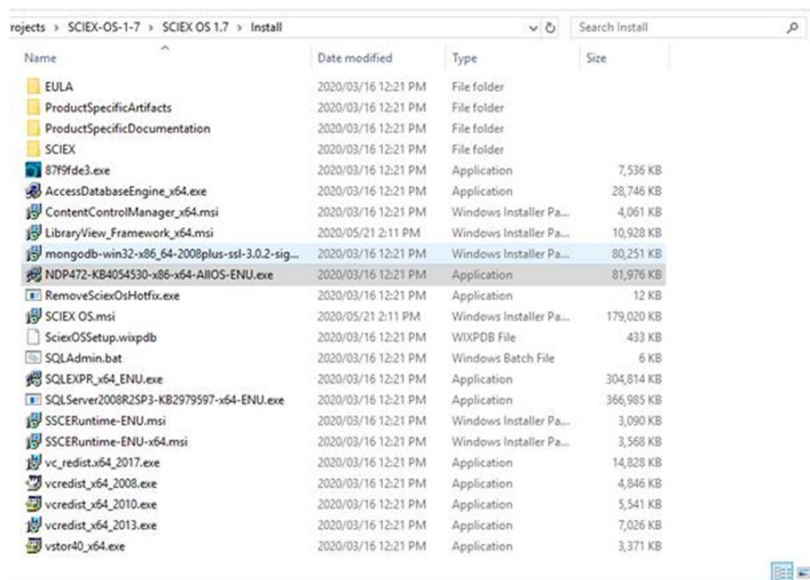

Release Notes

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate those manufacturers' products as supplied by SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products and are the sole and exclusive representations, warranties, and obligations of SCIEX. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser or for any adverse circumstances arising therefrom.

For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd., or their respective owners, in the United States and/or certain other countries. AB SCIEX™ is being used under license.

© 2020 DH Tech. Dev. Pte. Ltd.



AB Sciex Pte. Ltd.
Blk 33, #04-06 Marsiling Ind Estate Road 3
Woodlands Central Indus. Estate SINGAPORE 739256