

SCIEX OS 2.2 HotFix 1 Rev.2 Release Notes



Introduction

For information about the SCIEX OS 2.2 software, including requirements, supported firmware and configuration table versions, and known issues, refer to the *Release Notes* for SCIEX OS 2.2.

For information about previous software versions and HotFixes, refer to the *Release Notes* that came with the software version.

Where to Get Help

- *SCIEX OS 2.2 Software Release Notes*
- *SCIEX OS 2.2 Software Installation Guide*

New in SCIEX OS 2.2 HotFix 1 Rev.2

Fixed Issues

- The Installation Confirmation Tool failed after installation of SCIEX OS 2.2 HotFix 1. (BLT-3065)
- SCIEX OS 2.2 was not able to acquire data to the specified network folder using the Secure Network Account (SNA). (BLT-3526)
- When users tried to change the root directory to a network directory in SCIEX OS, the error message was shown: `The user does not have permission to create a folder.` (BLT-3486)
- A file read error was shown when users processed data with SCIEX OS while the Analyst software was acquiring data to the same file with the checksum on. (BLT-3578)

HotFixes and Patches Included

- N/A

Installation

Install the HotFix

Prerequisites
<ul style="list-style-type: none">• SCIEX OS 2.2 is installed. <hr/> <p>Note: SCIEX OS 2.2 HotFix 1 Rev.2 is compatible with SCIEX OS 2.2 HotFix 1 and can be installed on computers with HotFix 1 installed. HotFix 1 is removed silently during installation of this version.</p> <hr/>

1. Log on to the computer as a user with Administrator privileges.
2. Stop any acquisitions that are in progress and then deactivate the devices.
3. Close SCIEX OS.
4. Download **SCIEX OS 2.2 HotFix 1 Rev.2** from sciex.com/software-support/software-downloads.

Tip! To prevent potential installation issues, save the file to a location other than the computer desktop and then disconnect any external USB storage devices before starting the installation.

5. After the download is complete, right-click the `SCIEX-OS-2.2-HotFix-1-Rev.2.zip` file.
6. Click **Extract All**, browse to and select the destination folder, and then click **Extract**.
7. After the extraction is complete, browse to the extracted folder, and then double-click the `SCIEX_OS_2.2_HotFix_1_Rev.2.exe` file.
8. Follow the on-screen instructions to complete the installation.
9. Restart the computer.
10. Open SCIEX OS and then activate the devices.

Remove the HotFix

Remove the HotFix before removing SCIEX OS.

1. Log on to the computer as a user with Administrator privileges.
2. Deactivate the devices and then close SCIEX OS.
3. Close SCIEX OS.
4. Open Control Panel and then click **Programs and Features**.

5. Select **SCIEX OS 2.2 HotFix 1 Rev.2** and then click **Uninstall**.
6. Follow the on-screen instructions.

The HotFix is removed from the program list, and any files that were changed by the HotFix are reverted to the original SCIEX OS 2.2 software version.

7. Close Control Panel.

Updated Files

The SCIEX OS 2.2 HotFix 1 Rev.2 makes changes to the following folders and files.

C:\Program Files\SCIEX\SCIEX OS

- Clearcore.hsh
- Clearcore2.Data.Acquisition.Contracts.dll
- Clearcore2.SampleData.TransferQueue.dll
- OFX.Core.Contracts.dll
- OFX.Security.Library.dll
- ProductSpecificArtifacts.hsh
- Sciex.Clearcore.FMan.dll

C:\Program Files\SCIEX\SCIEX OS\Analytics\bin

- Analyst.NET.dll (unchanged, added to this location)
- Sciex.Clearcore.FMan.dll
- Sciex.MultiQuant.WiffFileState.exe

Known Issues and Limitations

- N/A

Contact Us

Customer Training

- In North America: NA.CustomerTraining@sciex.com
- In Europe: Europe.CustomerTraining@sciex.com
- Outside the EU and North America, visit sciex.com/education for contact information.

Online Learning Center

- [SCIEX Now Learning Hub](#)

SCIEX Support

SCIEX and its representatives maintain a staff of fully-trained service and technical specialists located throughout the world. They can answer questions about the system or any technical issues that might arise. For more information, visit the SCIEX website at sciex.com or contact us in one of the following ways:

- sciex.com/contact-us
- sciex.com/request-support

CyberSecurity

For the latest guidance on cybersecurity for SCIEX products, visit sciex.com/productsecurity.

Documentation

This version of the document supercedes all previous versions of this document.

To view this document electronically, Adobe Acrobat Reader is required. To download the latest version, go to <https://get.adobe.com/reader>.

To find software product documentation, refer to the release notes or software installation guide that comes with the software.

To find hardware product documentation, refer to the documentation DVD for the system or component.

The latest versions of the documentation are available on the SCIEX website, at sciex.com/customer-documents.

Note: To request a free, printed version of this document, contact sciex.com/contact-us.

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate those manufacturers' products as supplied by SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products and are the sole and exclusive representations, warranties, and obligations of SCIEX. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser or for any adverse circumstances arising therefrom.

(GEN-IDV-09-10816-D)

For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd., or their respective owners, in the United States and/or certain other countries (see sciex.com/trademarks).

AB Sciex™ is being used under license.

© 2022 DH Tech. Dev. Pte. Ltd.



AB Sciex Pte. Ltd.
Blk33, #04-06 Marsiling Industrial Estate Road 3
Woodlands Central Industrial Estate, Singapore 739256